U.S. English

*Product Number:
RQAFM1HS*

Reveal[®] Q+ HS

*for Aflatoxin M₁
(21–150 parts per trillion)*

For use with NEOGEN[®]'s Raptor[®] Integrated Analysis Platform

Refrigerate at 2–8°C (35–46°F) Do not freeze.

Reveal[®] Q+ HS for Aflatoxin M₁

Product Number: RQAFM1HS

The Toxin

Aflatoxin is a toxic and carcinogenic substance produced by certain strains of the molds *Aspergillus flavus* and *A. parasiticus*. When cows consume grain containing aflatoxin, they create a metabolite that is found in their milk, called aflatoxin M1.

Ingesting excessive amounts of the toxin can cause chronic health and performance problems or death in animals. Aflatoxin has been shown to cause liver damage or cancer, decreased milk, and egg production, immune system suppression, and interference with reproductive efficiency.

Intended Use and User

Reveal[®] Q+ HS for Aflatoxin M1 is intended for the screening of raw commingled bovine milk for aflatoxin M1. The test kit is designed for use by quality control personnel and others familiar with handling milk possibly contaminated with aflatoxin M1.

Assay Principles

Reveal Q+ HS for Aflatoxin M1 is a single-step lateral flow immunoassay based on a competitive immunoassay format. The milk is wicked through a reagent zone, which contains antibodies and receptors conjugated to colloidal gold particles. If aflatoxin M1 is present, it will be captured by the particle-antibody complex. The toxin-antibody-particle complex is then wicked onto a membrane that contains a test line. This test line captures any antibody gold particle complexes (in absence of aflatoxin M1) to concentrate and form a visible line. As the level of aflatoxin M1 in the sample increases, any free toxin will complex with the antibody gold particles. This allows less antibody-gold to be captured into the test line. Therefore, as the concentration of aflatoxin M1 in the sample increases, the test line density decreases. The membrane also contains a control line where an immune complex present in the reagent zone is captured by an antibody, forming a visible line. The control line will always form regardless of the presence of aflatoxin M1, ensuring the strip is functioning properly. This assay is meant to be performed using the Raptor[®] Integrated Analysis Platform.

Storage Requirements

Store kit components at temperatures between 2–8°C to ensure full shelf life. Test strips should remain capped in their original sample tube until used to ensure optimal performance.

Materials Provided

1. 1 stay-dry sample tube with 25 test strips
2. 25 disposable polypipettes

Materials Required, But Not Provided:

1. Raptor Integrated Analysis Platform (NEOGEN[®] item 9680)
2. Raptor cartridges (NEOGEN item 9681)

Precautions

1. The test strips must remain inside their stay-dry sample tube before use.
2. Store test kit at temperatures between 2–8°C when not in use. Do not freeze.
3. Do not use kit components beyond expiration date.

Test Preparation

1. Insert the test strip into the Raptor cartridge in any of the available slots.
2. Insert the cartridge containing the test strip into any of the 3 ports within the Raptor Integrated Analysis Platform.
 - a. The bar code on the test device will be read. If the QR code for the lot of devices has not been entered into the Raptor, the bar code reader in the front of the Raptor will turn on automatically. Scan the QR code found on the container storing the test devices.
 - b. The assay temperature for the Reveal Q+ HS for Aflatoxin M1 test is 65°C. This temperature is programmed in the Raptor.
 - c. When the cartridge is inserted into one of the 3 ports, the port will automatically begin to adjust to the proper temperature.
 - d. The user will not be able to proceed until the incubator temperature reaches 65°C.
3. Enter the sample ID by either scanning or using the on-screen keyboard.
4. Mix the milk sample prior to adding the sample into the cartridge.
5. Pipette 0.4 mL milk sample into the sample port located in the back of the cartridge. Press next. The analysis will take place automatically.
6. Insert the used pipette into the sample port located in the back of the cartridge. This will prevent double loading the same sample or loading a second sample into the same cartridge.
7. After a 5 minute incubation, the results will appear on the Raptor screen.

Interpretation of Results

The reader will use a lot-specific curve to calculate the results in parts per trillion (ppt) based on the line density values on the strip. Results reported less than 32 ppt are considered a negative result and results reported greater than 32 ppt and less than 53 ppt are a suspected positive result for the 50 ppt regulatory level and should be confirmed by alternative method such as Veratox for Aflatoxin M1. The range of quantification of the product is 53–150 ppt, therefore; results are fully quantitative from 53 ppt.

Performance Characteristics

Limit of detection: 21 ppt

Range of quantitation: 53–150 ppt

Customer Service

NEOGEN Customer and Technical Services can be contacted through [NEOGEN.com](https://www.NEOGEN.com) and product training is available by request.

SDS Information Available

Safety data sheets are available for all test kits at [NEOGEN.com](https://www.NEOGEN.com) or by calling 800.234.5333 or 517.372.9200.

Terms and Conditions

NEOGEN's full terms and conditions are available [online](#).

Warranty

NEOGEN makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, NEOGEN will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. NEOGEN shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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