

Reveal Procedure for Ruminant Feed

Note: Please read kit instructions completely before performing test.
Questions? Call 800/234-5333 or 517/372-9200.



1. Weigh 10 grams of sample and place in a Mason-type jar.



2. Add one level scoop of extraction additive.



3. Add 100 mL of extraction solution to the jar.



4. Heat in boiling water for 10 minutes.



5. Transfer 0.5 mL of extract to a sample tube.



6. Place a strip into the sample tube.



7. Interpret results after 15 minutes. If two lines appear, the sample is positive. If one line appears, the sample is negative.

Product Specifications

Testing time: Extraction boiling step: 10 minutes;
Test strip development: 15 minutes

Sensitivity: 1% in feed and feed supplements

Storage: Room temperature, 18-30°C



620 Leshar Place • Lansing, MI 48912
800/234-5333 (USA and Canada only) • 517/372-9200 • fax: 517/372-2006
e-mail: foodsafety@neogen.com • www.neogen.com