


Alert Procedure for Total Milk Allergen

Note: Please read kit instructions completely before performing test.
Questions? Call 800/234-5333 or 517/372-9200.

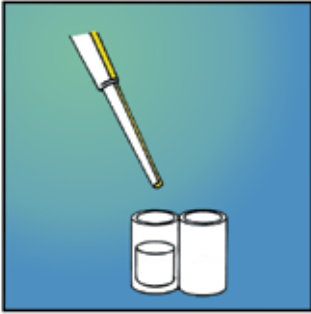

1. Choose one of the controls provided. Add 100 μ l from that control bottle to the first well.



2. Add 100 μ l or 3 drops of each sample extract to respective test well. Mix. Incubate for 10 minutes.



3. Dump liquid from wells.



4. Wash wells thoroughly 5 times with washing solution. Tap out liquid on absorbent paper towel.



5. Add 100 μ l of conjugate to each well. Mix. Incubate for 10 minutes.




6. Repeat steps 3-4 by dumping out the liquid, thoroughly washing the wells and tapping dry.



7. Add 3 drops of substrate to each well. Mix. Incubate for 10 minutes.



8. Add 3 drops of Red Stop to each well. Mix.

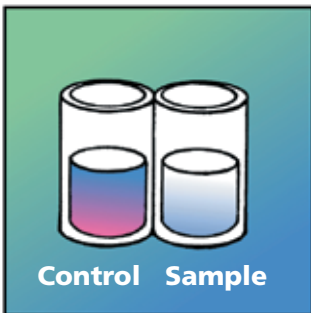

9. Visually read results. (Sample color lighter than the control indicates negative results.)

Product Specifications

- Controls provided: 5 ppm and 10 ppm
- Testing time: 30 minutes



620 Leshar Place • Lansing, MI 48912
800/234-5333 (USA and Canada only) • 517/372-9200 • fax: 517/372-2006
e-mail: foodsafety@neogen.com • www.neogen.com