*Read instructions carefully before using*

## for Histamine Extraction Kit

*Designed for use with Neogen's Veratox for Histamine test kit*

### INTENDED USE

Veratox for Histamine Extraction kit is intended for the quantitative analysis of histamine in scombroid species of fish, such as tuna, bluefish and mahi-mahi, fish meal, and dry animal protein.

### MATERIALS PROVIDED

1. 38 disposable extraction bottles (see **Precaution 1**)
2. 38 disposable filter syringes
3. 38 sample collection tubes, plastic (12 x 75 mm)
4. Directions for use

### MATERIALS REQUIRED BUT NOT PROVIDED

1. Blender
2. Graduated cylinder capable of measuring 90 mL (Neogen item 9368)
3. Timer (Neogen item 9426)
4. Test tube rack (Neogen item 9440)
5. Distilled or deionized water

### STORAGE

This kit may be stored at room temperature, 18–30°C (64–86°F).

### PRECAUTIONS

1. If using extraction bottles other than those included in the Veratox for Histamine Extraction Kit, make sure the bottles are plastic, **NOT GLASS**. Histamine adheres to glass, thus the use of glassware can affect test results.
2. Good laboratory practices should always be followed.
3. Always use a new extraction bottle, sample collection tube and filtering syringe for each sample to avoid cross-contamination.
4. Destroy and dispose of filtering syringe in a manner to prohibit reuse.
5. Do not take filtering material out of syringe. Material can cause skin and eye irritation.

### SAMPLE PREPARATION FOR EXTRACTION

#### 1. **Canned tuna: AOAC 937.07b**

Place entire contents of can (meat and liquid) into a blender. Blend until homogenous.

#### 2. **Fresh or thawed frozen raw fish: AOAC 937.07a**

Clean and eviscerate three fish. Cut three cross-sectional pieces 2.5 cm (1 inch) thick, from back of the pectoral fin, halfway to vent and one posterior to the vent. Debone slices and blend or grind combined samples until homogenous.

#### 3. **Dry Samples (example: Fishmeal)**

The sample to be tested should be collected according to accepted sampling techniques and be representative of the lot. After homogenization, grind part of the sample (minimum of 200 g) so at least 95% of the ground material passes through a 20 mesh sieve (about the particle size of fine espresso).

## SAMPLE EXTRACTION

**NOTE:** Glassware should not be used for extraction purposes as histamine may adhere to glass.

### Liquid/Wet Samples

1. Add **10 g** of the homogenous mixture to a clean disposable extraction bottle containing **90 mL** of distilled or deionized water.
2. Tightly cap and vigorously shake the bottle for **15-20 seconds** in order to suspend the fish tissue in the water.
3. Wait approximately **5 minutes**, then shake the bottle for **15-20 seconds** in order to re-suspend the fish tissue.
4. Wait an additional **5 minutes**, and again shake the bottle for **15-20 seconds** in order to re-suspend the fish tissue. Allow the tissue to settle to the bottom of the bottle for about **30 seconds**.
5. If necessary, filter the contents through folded filter paper or a histamine filter syringe (Neogen #9420H) into a clean container. The sample is now ready for testing with the Veratox for Histamine test kit. Proceed to the test kit's **SAMPLE EXTRACT DILUTION** step.

**ALTERNATIVE:** Centrifuge the sample and use the clear supernatant as the sample for extract dilution.

### Dry Samples

1. Add **10 g** of the homogenous mixture to a clean disposable extraction bottle containing **100 mL** of distilled or deionized water.
2. Tightly cap and vigorously shake the bottle for **15-20 seconds** in order to suspend the sample in the water.
3. Wait approximately **5 minutes**, then shake the bottle for **15-20 seconds** in order to re-suspend the sample.
4. Wait an additional **5 minutes**, and again shake the bottle for **15-20 seconds** in order to re-suspend the sample. Allow the sample to settle to the bottom of the bottle for about **30 seconds**.
5. If necessary, filter the contents through folded filter paper or a histamine filter syringe (Neogen #9420H) into a clean container. The sample is now ready for testing with the Veratox for Histamine test kit. Proceed to the test kit's **SAMPLE EXTRACT DILUTION** step.

**ALTERNATIVE:** Centrifuge the sample and use the clear supernatant as the sample for extract dilution.

## CUSTOMER SERVICE

Neogen Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all Neogen test kits, is available.

## MSDS INFORMATION AVAILABLE

Material safety data sheets (MSDS) are available for this test kit, and all of Neogen's test kits, on Neogen's website at [www.neogen.com](http://www.neogen.com), or by calling Neogen at 800/234-5333 or 517/372-9200.

## TERMS AND CONDITIONS

For Neogen's full terms and conditions, please visit [www.neogen.com/Corporate/termsconditions.html](http://www.neogen.com/Corporate/termsconditions.html).

## WARRANTY

Neogen Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, Neogen will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. Neogen shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

### North America Neogen Headquarters

620 Leshler Place, Lansing, MI 48912 USA  
800/234-5333 (USA/Canada) or 517/372-9200  
Fax: 517/372-2006 • [foodsafety@neogen.com](mailto:foodsafety@neogen.com)  
[www.neogen.com](http://www.neogen.com)

### Europe, Middle East and Africa Neogen Europe

The Dairy School, Auchincruive, Ayr  
KA6 5HW Scotland, UK  
+ 44 (0) 1292 525 600  
Fax: + 44 (0) 1292 525 601  
[info\\_uk@neogeneurope.com](mailto:info_uk@neogeneurope.com)  
[www.neogeneurope.com](http://www.neogeneurope.com)

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