Read instructions carefully before starting test

Reveal Q+

for T-2/HT-2

**For use with Neogen's line of AccuScan® readers,
including AccuScan Pro and AccuScan Gold**

THE TOXIN

T-2/HT-2 toxins are trichothecene mycotoxins produced by several species of *Fusarium* molds. Since T-2 toxin is readily metabolized to HT-2 toxin and the toxins have been shown to produce numerous adverse effects on many animals, these two mycotoxins are frequently evaluated together.

Animals affected by the toxins include swine, dairy cattle, poultry, dogs, cats and horses. Effects of the toxins include digestive disorders, hemorrhage, edema, oral lesions, dermatitis, and blood disorders. Damage caused by the toxins to the digestive track is irreversible. In the most severe cases, these toxins will cause death. T-2 toxin is the principal causal toxin in the human disease alimentary toxic aleukia.

Poultry studies have shown T-2 intoxication has led to a reduction in weight gain and other problems such as beak lesions, poor feathering, motor function impairment and increased susceptibility to *Salmonella* spp.

The best protection against these mycotoxins is monitoring for their presence in feeds and foods. That means testing all along the pathway from initial harvest of grains to the finished product.

INTENDED USE/USER

Reveal® Q+ for T-2/HT-2 is intended for the quantitative analysis of corn and corn products for T-2/HT-2 toxins. The test kit is designed for use by quality control personnel and others familiar with commodities possibly contaminated with T-2/HT-2 toxins.

ASSAY PRINCIPLES

Reveal Q+ for T-2/HT-2 is a single-step lateral flow immunochromatographic assay based on a competitive immunoassay format. The extract is wicked through a reagent zone, which contains antibodies specific for T-2/HT-2 toxins conjugated to colloidal gold particles. If T-2/HT-2 toxins are present, they will be captured by the particle-antibody complex. The T-2/HT-2 toxins-antibody-particle complex then is wicked onto a membrane, which contains a zone of T-2/HT-2 toxins conjugated to a protein carrier. This zone captures any uncomplexed T-2/HT-2 toxins antibody, allowing the particles to concentrate and form a visible line. As the level of T-2/HT-2 toxins in a sample increases, free T-2/HT-2 toxins will complex with the antibody-gold particles. This allows less antibody-gold to be captured in the test zone. Therefore, as the concentration of T-2/HT-2 toxins in the sample increases, the test line density decreases. Algorithms programmed into the reader convert these line densities into a quantitative result displayed in parts per billion (ppb). The membrane also contains a control zone where an immune complex present in the reagent zone is captured by an antibody, forming a visible line. The control line will always form regardless of the presence of T-2/HT-2 toxins, ensuring the strip is functioning properly.

STORAGE REQUIREMENTS

Store kit components at room temperature (18–30°C, 64–86°F) to ensure full shelf life. Test strips should remain capped in their original tubes until used to ensure optimal performance.

MATERIALS PROVIDED

1. 25 Reveal Q+ for T-2/HT-2 test strips
2. 25 red sample dilution cups
3. 25 clear sample cups
4. 1 bottle of sample diluent
5. Instructions for use

MATERIALS RECOMMENDED BUT NOT PROVIDED

1. Sample collection cups with lids (NEOGEN item 9428)
2. Agri-Grind grinder or equivalent (NEOGEN item 9401, 9453)
3. Scale capable of weighing 5–50 g ± 0.1 g (NEOGEN item 9427)
4. Timer (NEOGEN item 9426)
5. Reveal sample cup rack (NEOGEN item 9475)
6. AccuScan reader: AccuScan Pro or AccuScan Gold
7. Dispensing pump or graduated cylinder (NEOGEN item 9448, 9447)

8. Filter syringe (NEOGEN item 9420)
or
Centrifuge, mini (NEOGEN item 9330)
Microcentrifuge tubes (NEOGEN item 9172)
9. Sample collection tubes with caps (NEOGEN item 9421, 9421B)
10. Pipettor, 100 μ L (NEOGEN item 9272, 9278)
11. Pipette tips, 100 μ L (Neogen item 9407, 9410, 9417)
12. Pipettor, 500 μ L (Neogen item 9336)
13. Pipette tips, 200–1000 μ L (Neogen item 9464, 9487)

PRECAUTIONS

1. The test strips must remain inside the stay-dry tube before use.
2. Store test kit at room temperature (18–30°C, 64–86°F) when not in use.
Do not freeze.
3. Do not use kit components beyond expiration date.
4. Treat all used liquids, including sample extract, and labware as if contaminated with T-2/HT-2 toxins. Gloves and other protective apparel should be worn at all times.
5. To avoid cross-contamination, use clean glassware for each sample, and thoroughly wash all glassware between samples.
6. Ensure the device lot number and the curve details match the lot ID number selected on the reader. Failure to update the lot-specific QR code within the AccuScan Pro and AccuScan Gold reader will cause inaccurate results.

ACCUSCAN READER SET UP

AccuScan Pro reader

1. Enter the lot-specific QR code by selecting the QR code icon on the reader. Place the QR code into the cartridge and insert the cartridge into the reader.
NOTE: For instructions on manually entering sample IDs, see the AccuScan Pro user manual.
2. Return to the home screen and select the test strip icon. Touch the mycotoxin category, then select the **T-2/HT-2 Q+** test type.

AccuScan Gold reader

1. Enter the lot specific QR code by selecting **Scan QR** from the main screen. Place the lot specific QR code into the white cartridge adapter labeled Cal/QR and place the cartridge into the reader.
2. The valid code will be scanned by the reader and provide information on the lot number and expiry date. Verify this information is correct and then add the lot ID to the reader by pressing **Add Lot ID**.
NOTE: The lot ID for the current lot will now be stored with the test ID (e.g., T-2/HT-2) and can be selected when running a test.

PERFORMANCE CHARACTERISTICS

1. Limit of detection: 50 ppb
2. Range of quantitation: 50–600 ppb

NOTE: Samples greater than 600 ppb must be diluted and retested. Results below the range of quantitation should be reported as less than 50 ppb.

VALIDATED/VERIFIED MATRICES

Neogen continues to validate new commodities. Please contact a representative for the latest validated commodity list.

CUSTOMER SERVICE

NEOGEN Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all NEOGEN test kits, is available.

SDS INFORMATION AVAILABLE

Safety data sheets (SDS) are available for this test kit on NEOGEN's website at NEOGEN.com, or by calling NEOGEN at 800.234.5333 or 517.372.9200.

TERMS AND CONDITIONS

For NEOGEN's full terms and conditions, please visit www.NEOGEN.com/en/terms-and-conditions.

WARRANTY

NEOGEN Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, NEOGEN will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. NEOGEN shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

TEST KITS AVAILABLE FROM NEOGEN

Natural toxins

- Aflatoxin, DON, ochratoxin, zearalenone, T-2/HT-2 toxins, fumonisin, histamine

Foodborne bacteria

- *E. coli* O157:H7, *Salmonella*, *Listeria*, *Listeria monocytogenes*, *Campylobacter*, *Staphylococcus aureus*, *Salmonella enteritidis*

Sanitation

- ATP, yeast and mold, total plate count, generic *E. coli* and total coliforms, protein residues

Food allergens

- Almonds, crustaceans, eggs, gliadin, hazelnut, milk, mustard, peanuts, sesame, soy, walnuts, multi-treenut

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