Read instructions carefully before starting test

Reveal[®] Q+ MAX
for Zearalenone
Quantitative Test
With Aqueous Extraction

THE TOXIN

Zearalenone is primarily produced by the mold *Fusarium graminearum*, which also commonly produces deoxynivalenol (DON). Hence, there is evidence that if zearalenone is detected, there is a high probability that other fusarial mycotoxins may be present. Zearalenone is classified as an estrogenic mycotoxin because it frequently causes estrogenic responses in animals.

When zearalenone-contaminated feed or grain is eaten by livestock, it can cause a wide variety of reproductive problems. In swine, it causes vulvovaginitis, low birth weights, fetal reabsorption, aborted pregnancies, reduced litter sizes, abnormal estrus and feminization of immature males. Zearalenone can delay the breeding process and cost the producer significant economic and physical losses.

Livestock producers are becoming increasingly aware of zearalenone problems and have looked for ways to reduce risks related to contaminated feed.

The best protection against mycotoxins is monitoring for their presence in feeds and foods. That means testing all along the pathway from initial harvest of grains to the finished product.

The FDA has not issued any advisory levels for zearalenone.

European Union regulations for zearalenone are as follows:

Foodstuffs	Maximum levels parts per billion (ppb)
Unprocessed corn	200
Other unprocessed cereals and corn for direct human consumption	100
Other cereals for direct human consumption	75
Bread, pastries, breakfast cereals, snacks, corn-based snacks and corn-based breakfast cereals	50
Baby foods and cereals intended for infants	20

INTENDED USE/USER

Reveal® Q+ MAX for Zearalenone is intended for the quantitative analysis of corn, wheat and other commodities for zearalenone. This test utilizes an aqueous extraction, thus eliminating the need for the disposal of hazardous waste. The test kit is designed for use by quality control personnel and others familiar with commodities possibly contaminated by zearalenone.

ASSAY PRINCIPLES

Reveal Q+ MAX for Zearalenone is a single-step lateral flow immunoassay. In the test, the extract is wicked through the reagent zone, which contains antibodies specific for zearalenone conjugated to colloidal gold particles (gold complex). If zearalenone is present, it will be captured by the gold complex. The gold complex, along with any free gold-complex, is then wicked onto a membrane, which contains a zone of zearalenone conjugated to a protein carrier. This zone captures any unbound gold complex, allowing the particles to concentrate and form a visible line.

As the level of zearalenone in a sample increases, free zearalenone will bind with the gold complex allowing less gold complex to be captured in the test zones. Therefore, as the concentration of zearalenone in the sample increases, the test line density decreases. Algorithms programmed into the AccuScan® readers convert the line intensities into a quantitative result displayed in parts per billion (ppb). The membrane also contains a control zone where an immune complex present in the reagent zone is captured by an antibody, forming a visible line. The control line will always form regardless of the presence of zearalenone, ensuring the strip is functioning properly.

STORAGE REQUIREMENTS

Store kit components at room temperature between 18–30°C (64–86°F) to ensure full shelf life. Test strips should remain capped in their original tubes until used to ensure optimal performance.

MATERIALS PROVIDED

1. 25 Reveal Q+ MAX for Zearalenone test strips
2. 25 red conical sample dilution cups
3. 25 clear sample cups
4. 2 bottles of red-labeled sample diluent
5. 1 bottle of yellow-labeled dilution buffer (for samples > 800 ppb)
6. 25 MAX 1 Aqueous Extraction packets
7. Instructions for use

MATERIALS RECOMMENDED BUT NOT PROVIDED

1. Extraction Materials
 - a. Sample collection cups with lids 125 mL (Neogen item 9428, 9428B)
 - b. Sample collection tubes with caps (Neogen item 9421, 9421B)
 - c. Filter syringes, Whatman #4 filter paper or equivalent (Neogen item 9420, 9429, 9519)
 - d. Dispensing pump or graduated cylinder (Neogen item 9448, 9447)
2. Agri-Grind grinder or equivalent (Neogen item 940, 9453)
3. Scale capable of weighing 5–50 grams \pm 0.1 g (Neogen item 9427)
4. Timer (Neogen item 9426, 9452)
5. Reveal sample cup rack (Neogen item 9475)
6. Pipettor, 100 μ L (Neogen item 9860, 9272)
7. Pipettor 100–1000 μ L (Neogen item 9463)
8. Pipettor, 250 μ L (Neogen item 9692)
9. Pipette tips, 100 μ L–1 mL (Neogen item 9464, 9487)
10. Pipette tips, 1–200 μ L (Neogen item 9407, 9410, 9417)
11. Distilled or deionized water
12. AccuScan Gold reader (Neogen item 9595), AccuScan Pro reader (Neogen item 9565) or Raptor® Integrated Analysis Platform (Neogen item 9680)
13. Raptor cartridges (Neogen item 9681)
14. Raptor 400 μ L Exact Volume pipettes (Neogen item 9682)

PRECAUTIONS

1. The test strips must remain inside the stay-dry tube before use.
2. Store test kit at room temperature 18–30°C (64–86°F) when not in use. Do not freeze.
3. Do not use kit components beyond expiration date.
4. Treat all used liquids, including sample extract, and labware as if contaminated with zearalenone.
5. Ensure the device lot number and the curve details match the lot ID number selected on the reader. Failure to update the lot-specific QR code within the AccuScan reader will cause inaccurate results.
6. Commodity extracts should have a pH of 6–8 before testing. Excessively acidic or alkaline samples should be adjusted. For instructions on adjusting pH contact a Neogen representative or Technical Services.

ACCUSCAN GOLD READER SETUP

1. Enter the lot-specific QR code by selecting **Scan QR** from the main screen. Place the lot specific QR code into the white cartridge adapter labeled Cal/QR and place the cartridge into the reader.
2. The valid code will be scanned by the reader and provide information on the lot number and expiry date. Verify this information is correct and then add the lot ID to the reader by pressing **Add Lot ID**.

NOTE: The lot ID for the current lot will now be stored with the test ID (e.g., zearalenone) and can be selected when running a test.

ACCUSCAN PRO READER SETUP

Enter the lot-specific QR code by selecting the QR code icon on the reader. Place the QR code into the cartridge and insert the cartridge into the reader.

NOTE: For instructions on manually entering sample IDs, see the AccuScan Pro user manual.

SAMPLE PREPARATION

The sample to be tested should be collected according to accepted sampling techniques (see FGIS sampling protocol or contact your Neogen representative). Obtain a representative sample (minimum 100 g). Grind the sample so at least 95% of the ground material passes through a 20 mesh sieve (about the particle size of fine espresso).

TESTING PROTOCOL FOR WHEAT

Sample Extraction

1. Weigh out 10 ± 0.1 g of sample into extraction cup.
2. Add contents of one (1) MAX 1 aqueous extraction packet to the extraction cup.
3. Add **50 mL** distilled or deionized water to the extraction cup.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman No. 4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

SAMPLE Extraction (FGIS Method)

1. Weigh out 50 ± 0.1 g of sample into extraction container.
2. Add contents of one (1) MAX 1-G50 aqueous extraction packet to the extraction cup.
3. Add **250 mL** distilled or deionized water to the extraction container.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman #4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

Testing Procedure

1. Place the appropriate number of red sample dilution cups and clear sample cups into a sample cup rack. Label cups if necessary.
2. Add 100 μ L of sample diluent (red-labeled bottle) to each red sample dilution cup.
3. Add 100 μ L of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
4. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
5. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
6. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5 minute incubation.
7. **For the AccuScan Gold reader:** **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: Q+M ZEAR Curve 1
For the AccuScan Pro reader: **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: ZEARQ+MAX Curve 1

Dilution Procedure

Samples **greater than 500 ppb** will need to be diluted and re-tested.

1. Add 100 μ L sample filtrate to sample collection tube.
2. Add 200 μ L dilution buffer (yellow-labeled bottle) to the sample collection tube. Mix by pipetting up and down 5 times.
3. Add 100 μ L of sample diluent (red-labeled bottle) to a red sample dilution cup.
4. Add 100 μ L of diluted sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
5. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
6. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
7. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5-minute incubation
8. **For the AccuScan Gold reader:** **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: Q+M ZEAR Curve 1
For the AccuScan Pro reader: **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: ZEARQ+MAX Curve 1
9. Multiply your result by 3.

TESTING PROTOCOL FOR CORN

Sample Extraction

1. Weigh out 10 ± 0.1 g of sample into extraction cup.
2. Add contents of one (1) MAX 1 aqueous extraction packet to the extraction cup.
3. Add **50 mL** distilled or deionized water to the extraction cup.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman No. 4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

Sample Extraction (FGIS Method)

1. Weigh out 50 ± 0.1 g of sample into extraction container.
2. Add contents of one (1) MAX 1-G50 aqueous extraction packet to the extraction cup.
3. Add **250 mL** distilled or deionized water to the extraction container.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman #4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

Testing Procedure

1. Place the appropriate number of red sample dilution cups and clear sample cups into a sample cup rack. Label cups if necessary.
2. Add 100 μ L of sample diluent (red-labeled bottle) to each red sample dilution cup.
3. Add 100 μ L of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
4. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
5. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
6. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5 minute incubation.
7. **For the AccuScan Gold reader: Select CATEGORY: Mycotoxin Q+MAX
TEST NAME: Q+M ZEAR Curve 2**
**For the AccuScan Pro reader: Select CATEGORY: Mycotoxin Q+MAX
TEST NAME: ZearQ+MAX Curve 2**

Dilution Procedure

Samples **greater than 500 ppb** will need to be diluted and re-tested.

1. Add 100 μ L sample filtrate to sample collection tube.
2. Add 200 μ L dilution buffer (yellow-labeled bottle) to the sample collection tube. Mix by pipetting up and down 5 times.
3. Add 100 μ L of sample diluent (red-labeled bottle) to a red sample dilution cup.
4. Add 100 μ L of diluted sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
5. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
6. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
7. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5 minute incubation.
8. **For the AccuScan Gold reader: Select CATEGORY: Mycotoxin Q+MAX
TEST NAME: Q+M ZEAR Curve 2**
**For the AccuScan Pro reader: Select CATEGORY: Mycotoxin Q+MAX
TEST NAME: ZearQ+MAX Curve 2**
9. Multiply your result by 3.

TESTING PROTOCOL FOR CORN GLUTEN MEAL

Sample Extraction

1. Weigh out 10 ± 0.1 g of sample into extraction cup.
2. Add contents of one (1) MAX 1 aqueous extraction packet to the extraction cup.
3. Add **50 mL** distilled or deionized water to the extraction cup.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman No. 4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

Sample Extraction (FGIS Method)

1. Weigh out 50 ± 0.1 g of sample into extraction container.
2. Add contents of one (1) MAX 1-G50 aqueous extraction packet to the extraction cup.
3. Add **250 mL** distilled or deionized water to the extraction container.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman #4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

Testing Procedure

1. Place the appropriate number of red sample dilution cups and clear sample cups into a sample cup rack. Label cups if necessary.
2. Add 200 μ L of sample diluent (red-labeled bottle) to each red sample dilution cup.
3. Add 300 μ L of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
4. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
5. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
6. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5 minute incubation.
7. **For the AccuScan Gold reader:** **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: Q+M ZEAR Curve 2
For the AccuScan Pro reader: **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: ZearQ+MAX Curve 2

Dilution Procedure

Samples **greater than 500 ppb** will need to be diluted and re-tested.

1. Add 200 μ L sample filtrate to sample collection tube.
2. Add 400 μ L dilution buffer (yellow-labeled bottle) to the sample collection tube. Mix by pipetting up and down 5 times.
3. Add 200 μ L of sample diluent (red-labeled bottle) to a red sample dilution cup.
4. Add 300 μ L of diluted sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times.
5. Transfer 100 μ L of diluted sample extract into a new **clear** sample cup.
6. Place a new Reveal Q+ MAX for Zearalenone test strip with the sample end down into the clear sample cup and set timer for **5 minutes**. Ensure the test strip comes into contact with liquid and begins to wick.
7. Remove the strip from the sample cup after it has developed for **5 minutes**. Test strips must be read within 30 seconds of completion of the 5-minute incubation.
8. **For the AccuScan Gold reader:** **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: Q+M ZEAR Curve 2
For the AccuScan Pro reader: **Select CATEGORY: Mycotoxin Q+MAX**
TEST NAME: ZearQ+MAX Curve 2
9. Multiply your result by 3.

READING TEST RESULTS

NOTE: Test strips should be read within **30 seconds** of completion of the 5-minute incubation. Refer to **AccuScan Reader Set Up** for test selection and set up information.

1. Select the assay type (e.g., zearalenone) from the menu and ensure the device lot number matches the lot ID number selected on the reader.

NOTE: Failure to update the lot-specific QR code will cause inaccurate results.

2. Fully insert the Reveal Q+ test strip into the R-labeled cartridge adapter with the sample end first and results facing out.

3. Insert the cartridge with test strip upside-down into the AccuScan Gold reader (the test lines will face downward into the reader). If using the AccuScan Pro reader, insert the cartridge right side up. The reader will automatically begin analyzing the cartridge.

CAUTION: Removing cartridge prior to completion can result in invalid readings.

4. The AccuScan reader will analyze the test strip and results will be displayed and stored in the reader.


NOTES:

1. Ensure device is fully inserted into cartridge.
2. Readings should be made between **5 and 5.5 minutes**. Readings after 5.5 minutes may be inaccurate due to over-development of the device.

TEST PROCEDURE — RAPTOR INTEGRATED ANALYSIS PLATFORM

Sample Extraction

1. Weigh out 10 ± 0.1 g of sample into extraction cup.
2. Add contents of one (1) MAX 1 aqueous extraction packet to the extraction cup.
3. Add **50 mL** distilled or deionized water to the extraction cup.
4. Vigorously shake, using hand or mechanical means, for 3 minutes, or blend for 1 minute.
5. Allow the sample to settle, then filter with a filter syringe or Whatman No. 4 filter paper, collecting a minimum of 3 mL filtrate into a sample collection tube.

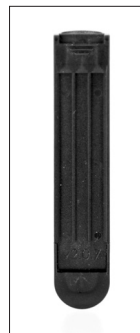Testing Procedure — Corn or wheat

1. Place the appropriate number of red sample dilution cups into a sample cup rack. Label cups if necessary.
2. Add 250 μ L of sample diluent (red-labeled bottle) to each red sample dilution cup.
3. Add 250 μ L of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times. Continue to step 4 below.

Testing Procedure — Corn gluten meal

1. Place the appropriate number of red sample dilution cups and clear sample cups into a sample cup rack. Label cups if necessary.
2. Add 200 μ L of sample diluent (red-labeled bottle) to each red sample dilution cup.
3. Add 300 μ L of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times. Continue to step 4 below.
4. Fully insert a Reveal Q+ MAX for Zearalenone test strip into a Raptor cartridge.
5. Insert the Raptor cartridge containing the test strip into any of the three ports within the Raptor.
 - a. The bar code on the test strip will be read. The system identifies the type of test strip and the lot number. If the lot number is not found in the system, the bar code reader on the front of the Raptor will turn on automatically.
 - b. Scan the QR code found on the tube containing the test strips. The information will be stored on the reader. (The QR code can also be found on the kit Certificate of Analysis.)
6. Enter the Sample ID if desired.
7. Select **Curve 1** for wheat or **Curve 2** for corn and corn gluten meal samples.
8. Add 400 μ L of sample from the red sample dilution cup to the Raptor cartridge.
 - a. The Raptor system will start automatically.
 - b. Additional samples can be started in the other ports while the first sample is processing.
9. Results will be displayed on the Raptor screen after the 5-minute incubation is complete.

DILUTION PROCEDURE — RAPTOR INTEGRATED ANALYSIS PLATFORM

Samples **greater than 500 ppb** must be diluted and re-tested.

Dilution Procedure — Corn or wheat

1. Add 200 μL of sample filtrate to a sample collection tube.
2. Add 400 μL of dilution buffer (yellow-labeled bottle) to the sample collection tube. Mix by pipetting up and down 5 times.
3. Add 250 μL of sample diluent (red-labeled bottle) to a red sample dilution cup.
4. Add 250 μL of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times. Continue to step 5 in the following section.

Dilution Procedure — Corn gluten meal

1. Add 200 μL of sample filtrate to a sample collection tube.
2. Add 400 μL dilution buffer (yellow-labeled bottle) to the sample collection tube. Mix by pipetting up and down 5 times.
3. Add 200 μL of sample diluent (red-labeled bottle) to a red sample dilution cup.
4. Add 300 μL of sample extract to the red sample dilution cup containing the sample diluent. Mix by pipetting up and down 5 times. Continue to step 5 below.
5. Fully insert a Reveal Q+ MAX for Zearalenone test strip into a Raptor cartridge.
6. Insert the Raptor cartridge containing the test strip into any of the three ports within the Raptor.
 - a. The bar code on the test strip will be read. The system identifies the type of test strip and the lot number. If the lot number is not found in the system, the bar code reader on the front of the Raptor will turn on automatically.
 - b. Scan the QR code found on the tube containing the test strips. The information will be stored on the reader. (The QR code can also be found on the kit Certificate of Analysis.)
7. Enter the Sample ID if desired.
8. Select **Curve 1** for wheat or **Curve 2** for corn and corn gluten meal samples.
9. Add 400 μL of sample from the red sample dilution cup to the Raptor cartridge.
10. The Raptor system will start automatically.
11. Additional samples can be started in the other ports while the first sample is processing.
12. Results will be displayed on the Raptor screen after the 5- minute incubation is complete.
13. Multiply your result by 3.

PERFORMANCE CHARACTERISTICS

Corn:

Limit of detection: 36 ppb

Range of detection: 50–500 ppb

Wheat:

Limit of detection: 21 ppb

Range of detection: 25–500 ppb

VALIDATED MATRICES

Wheat, corn and corn gluten meal.

NOTE: Neogen continues to validate new commodities. Please contact a representative for the latest validated commodity list.

CUSTOMER SERVICE

Neogen Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all Neogen test kits, is available.

SDS INFORMATION AVAILABLE

Safety data sheets (SDS) are available for this test kit, and all of Neogen's test kits, on Neogen's website at foodsafety.neogen.com, or by calling Neogen at 800/234-5333 or 517/372-9200.

TERMS AND CONDITIONS

For Neogen's full terms and conditions, please visit www.neogen.com/en/terms-and-conditions.

WARRANTY

Neogen Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, Neogen will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. Neogen shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

TESTING KITS AVAILABLE FROM NEOGEN

Natural toxins

- Aflatoxin, DON, ochratoxin, zearalenone, T-2/HT-2 toxins, fumonisin, histamine

Foodborne bacteria

- *E. coli* O157:H7, *Salmonella*, *Listeria*, *Listeria monocytogenes*, *Campylobacter*, *Staphylococcus aureus*, *Salmonella enteritidis*

Sanitation

- ATP, yeast and mold, total plate count, generic *E. coli* and total coliforms, protein residues

Food allergens

- Almonds, crustaceans, eggs, gliadin, hazelnut, milk, mustard, peanuts, sesame, soy, walnuts, multi-treenut

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