Read instructions carefully before starting test

Fluorescence Standards for ANSR Reader Certification

INTENDED USE

The ANSR[®] fluorescence standards provide for the certification of the performance of the ANSR reader.

MATERIALS PROVIDED

1. 16 reaction tubes of fluorescence standard
2. 16 reaction tubes of blank standard
3. 1 bottle of lysis reagent suspension buffer
4. Permanent caps
5. 1 kit insert
6. 2 foam-tipped swabs

EQUIPMENT REQUIRED

1. ANSR reader (Neogen item 9828)
2. Computer and software for connection to ANSR reader (Neogen item 9832)
3. Pipettor, 10-100 uL, 8-channel (Neogen item 9388)
4. Pipette tips, 100 uL, sterile, filtered (Neogen item 9389)
5. Vortex, adjustable (Neogen item 9494)
6. Pipettes, sterile serological 10 mL (Neogen item 8686)
7. Pipette pump for serological tubes, 10 mL (Neogen item 9277)

OTHER MATERIALS REQUIRED

1. Reagent reservoir
2. Strip tube rack
3. 70% isopropanol
4. Lint-free wipes

STORAGE

Store ANSR reagents at 2–8°C. After removing reaction tubes from the foil pouch, promptly re-seal the pouch. Leave the desiccant pack in the pouch at all times.

PRECAUTIONS

1. Use good microbiology laboratory practices.
2. **Dispose of used pipette tips in a covered container containing a fresh solution of 10% bleach. The 10% bleach solution should be made fresh each day. Use 9 parts water with 1 part commercially available household-strength bleach to make the 10% bleach solution. Undiluted stock solutions of bleach should be used within 30 days of opening.**
3. **Discard bleach solution and tips as regular waste at the end of each day.**
4. Do not use reagents beyond the expiration date.
5. Remove reaction tubes from the foil pouch just before use. Reseal the pouch to avoid prolonged exposure to light.
6. Do not remove caps from reaction tubes after the assay has been started in order to prevent accidental contamination of the environment with amplification products.
7. Exercise care in all pipetting steps to avoid cross-contamination of samples.
8. Complete all assay steps in sequence, avoiding delays between steps.
9. Tap reaction tubes on bench top to make sure lyophilized reagents are at the bottom of the tube prior to adding lysis reagent suspension buffer.

PRE-TEST CLEAN-UP

1. Put on a disposable set of gloves before handling the ANSR reader.
2. Clear space to clean the reader at least 3 meters away from open flames, as isopropanol is flammable.
3. Power down the reader and allow to cool.
4. Wipe all surfaces with lint-free wipes wetted with isopropanol. Use a sufficient amount of alcohol so that all surfaces are visibly wetted during the process.
5. Prepare a small aliquot of isopropanol.
6. Wet a foam-tipped swab (included) with the aliquot of isopropanol and gently clean the tube wells of the ANSR reader. Use a new swab for each set of 8 wells. Allow instrument to air-dry for 5 minutes.
7. Dispose of used materials and gloves.

ANSR TEST PROCEDURE

1. Remove the kit from the refrigerator and allow it to warm at room temperature (15–25°C) for 15 minutes.
2. Connect the ANSR reader to the computer via USB or Ethernet and turn the computer on.
3. Turn on the ANSR reader. The reader will preheat to $56 \pm 1^\circ\text{C}$.
4. Start the ANSR software and click the connect button. Confirm that the computer connects to the reader.
5. Click Run and select the appropriate test type from the drop-down menu (9847 ANSR Standards, 9847 Blank Standards).
6. Enter user information, kit lot number and sample identifiers (1–16).
NOTE: For instructions on using the reader and software, see the user guide that came with the ANSR reader.
7. Remove the appropriate reaction tube strips from the foil pouch and place in a reader.
8. Dispense approximately 3 mL of lysis reagent suspension buffer into a multi-channel pipette reagent reservoir.
9. Remove the caps from the reaction tube strips.
10. Using an 8-channel pipette and 100 μL filtered tips, transfer 50 μL of the lysis reagent suspension buffer to the reaction tubes. Place the provided permanent caps on the reaction tubes.
11. Vortex the reaction tubes at medium/high speed (about 2500 RPM) for 2–3 seconds and place them in the ANSR reader.
12. Close the lid and click Start in the ANSR software to begin the assay.
13. Repeat steps 5–12 for the 9847 Blank Standards, using the ANSR blank standard tubes.
NOTE: Fluorescence standards and blank standards need to be run and pass to certify the reader.

INTERPRETATION OF RESULTS

The ANSR software will first conduct a check of the control channel (HEX) to determine if the assay is valid. If the control is out of specification, an invalid will result for all channels.

If the control channel is determined to be valid, the ANSR software will then indicate the test results as positive or negative for the fluorescence performance of the test channels (FAM/ROX).

For the 9847 ANSR Standards assay:

If any negative or invalid results are produced, immediately perform the following:

1. Start a new assay in the ANSR software and re-select the 9847 test type.
2. Re-enter user information, kit lot number and sample identifiers (1–16).
3. Remove the tube strip from the reader, vortex for 2–3 seconds, then return to the reader.
4. Close the lid and click Start to repeat the assay.

Depending on the ANSR unit, and the last time the unit was synchronized to the ANSR Standards, it may take up to three attempts for the unit to pass. If the unit does not pass after three (3) attempts using the same reconstituted 9847 standards, the unit is not certified. Contact Neogen Technical Service for further information.

For the 9847 Blank Standards assay:

If any positive results are produced, immediately perform steps 1–4 listed above.

DISPOSAL

Do not remove permanent caps from the ANSR reaction tubes, even when disposing of them. Reaction tubes can be disposed of as non-biohazardous waste. It is recommended that they be placed in sealable plastic bags and immediately disposed of to protect against accidental opening.

CUSTOMER SERVICE

Neogen Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all Neogen test kits, is available.

SDS INFORMATION AVAILABLE

Safety data sheets (SDS) are available for this test kit, and all of Neogen's test kits, on Neogen's website at foodsafety.neogen.com, or by calling Neogen at 800/234-5333 or 517/372-9200.

TERMS AND CONDITIONS

Please visit <http://www.neogen.com/en/terms-and-conditions> for Neogen's full terms and conditions.

WARRANTY

Neogen Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, Neogen will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. Neogen shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

PRODUCTS AVAILABLE

- 9465 **Environmental sampling sponges** (with Dey Engley (DE) neutralizing broth) – pack of 100
- 36002 **Sampling sponges, pre-moistened, HiCap** – pack of 100
- 9822 **ANSR for *E. coli* O157:H7** – up to 96 tests
- 9824 **ANSR for *Listeria monocytogenes*** – up to 96 tests
- 9870 **ANSR for *Salmonella*** – up to 96 tests
- 9871 **ANSR for *Listeria*** – up to 96 tests
- 9872 **ANSR for *Campylobacter*** – up to 96 tests
- 9873 **ANSR for *Listeria Right Now*[™]** – up to 96 tests
- 9837 **Complete ANSR Equipment Kit** – Reader, computer, pipette set, dry bath heaters and inserts, thermometers, vortex, timer, rack

North America Neogen Headquarters

800/234-5333 (USA/Canada)
foodsafety@neogen.com
foodsafety.neogen.com

Europe, Middle East and Africa Neogen Europe

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China Neogen Bio-Scientific Technology

+86 21 6271 7013
info@neogenchina.com.cn
www.neogenchina.com.cn

India Neogen Food and Animal Security

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