

The logo for Soleris, featuring the word "soleris" in a bold, lowercase, sans-serif font. A stylized, curved line or swoosh passes through the letters, starting under the 's' and ending under the 's'.

Yeast and Mold Supplement Part #YMY-110Y

10 mL per vial – 4 vials per box

Each Vial Contains:

Chlortetracycline..... 100 mg

Usage Instructions:

1. For use with the DYM-109C Soleris vial
2. Add 10 mL of sterile water to the vial and gently mix
3. Add 0.3 mL of reconstituted supplement to each DYM-109C vial
4. Gently invert vial 3 times to thoroughly mix contents
5. The supplement should be used within 7 days following reconstitution. Store refrigerated.



620 Leshar Place, Lansing, MI 48912
800/234-5333 (USA/Canada) or 517/372-9200 • fax: 517/372-2006
e-mail: foodsafety@neogen.com • www.neogen.com