Read instructions carefully before starting test

Reveal[®]

for Aflatoxin Ethanol Extraction

Bulk Kit — 100 tests

THE TOXIN

Aflatoxin is a toxic and carcinogenic substance produced by certain strains of the molds *Aspergillus flavus* and *A. parasiticus*. There are four principle types of aflatoxin: B₁, B₂, G₁ and G₂. Aflatoxin B₁ is the most frequently encountered of the group and the most toxic. The commodities most affected by aflatoxin are corn, peanuts, cottonseed, milo, and the majority of tree nuts.

The affects in animals of ingesting excessive amounts of the toxin range from chronic health and performance problems to death. Aflatoxin has been shown to cause liver damage or cancer, decreased milk and egg production, immune suppression and interference with reproductive efficiency.

The FDA has set maximum allowable levels of aflatoxin in food and feed. Therefore, accurate determination of the presence of the toxin is of major importance to those monitoring the quality of food and feed in which aflatoxin may occur. Testing these commodities for the toxin requires careful sampling, chemical extraction, sanitation and quantitative analysis.

The FDA has issued regulatory levels for aflatoxin as follows:

For	Level	Commodities
Humans	20 ppb	All food except milk
All animal species	20 ppb	All feed (exceptions below)
Exceptions:		
Breeding cattle, breeding swine, mature poultry	100 ppb	Corn
Finishing swine (> 100 lbs.)	200 ppb	Corn
Finishing beef cattle	300 ppb	Corn
Finishing beef cattle, swine, poultry	300 ppb	Cottonseed meal

Reveal[®] for Aflatoxin is intended for the screening of corn for aflatoxin. The test kit is designed for use by quality control personnel and others familiar with commodities possibly contaminated by aflatoxin.

ASSAY PRINCIPLES

Reveal for Aflatoxin is a single-step lateral flow immunochromatographic assay based on a competitive immunoassay format. The extract is wicked through a reagent zone, which contains antibodies specific for aflatoxin conjugated to colloidal gold particles. If aflatoxin is present, it will be captured by the particle-antibody complex. The aflatoxin-antibody-particle complex then is wicked onto a membrane, which contains a zone of aflatoxin conjugated to a protein carrier. This zone captures any uncomplexed aflatoxin antibody, allowing the particles to concentrate and form a visible line. If aflatoxin is present above a pre-determined limit, no line will form. Therefore, a positive sample will result in the absence of a line in the test zone. Alternatively, a negative sample will form a visible line in the test zone. The membrane also contains a control zone where an immune complex present in the reagent zone is captured by an antibody, forming a visible line. The control line will always form regardless of the presence of aflatoxin, ensuring the strip is functioning properly.

STORAGE REQUIREMENTS

Store kit components at room temperature (18-30°C, 64-86°F) to assure full shelf life. Test strips should remain capped in their original sample tubes until used to assure optimal performance.

MATERIALS PROVIDED

1. 100 aflatoxin test strips
2. 100 sample cups
3. 4 bottles of sample diluent

MATERIALS RECOMMENDED BUT NOT PROVIDED

1. 65% ethanol solution
2. Distilled or deionized water
3. Sample collection cups with lids (Neogen item #9428)
4. Agri-Grind grinder or equivalent (Neogen item #9401)
5. Scale capable of weighing 10 grams \pm 0.1 g (Neogen item #9427)
6. Timer (Neogen item #9426)
7. Reveal sample cup rack (Neogen item #9475)
8. Cylinder, 50 mL graduated (Neogen item #9447)
9. Pipettor, 100 μ L (Neogen item 9272, 9860)
10. Pipette tips, 100 μ L (Neogen item 9407, 9410, 9417)
11. Pipettor, 500 μ L (Neogen item 9336, 9291)
12. Pipette tips, 200–1000 μ L (Neogen item 9464, 9487, 9293, 9292)
13. Reveal AccuScan Gold Reader

PRECAUTIONS

1. The test strips must remain inside the stay-dry tube before use.
2. Ethanol solution is highly flammable. Keep container tightly closed, and away from heat, sparks, open flame and those smoking. It is toxic if swallowed, or if vapor is inhaled. Avoid contact with skin.
3. Store test kit at room temperature (18-30°C, 64-86°F) when not in use, do not freeze.
4. Do not use kit components beyond expiration date.
5. Treat all used liquids, including sample extract, and labware as if contaminated with aflatoxin. Gloves and other protective apparel should be worn at all times.
6. To avoid cross-contamination, use clean glassware for each sample, and thoroughly wash all glassware between samples.

SAMPLE PREPARATION AND EXTRACTION

The sample to be tested should be collected according to accepted sampling techniques (see GIPSA sampling protocol or contact your Neogen representative). Obtain a representative sample (minimum 100 g). Grind the sample so at least 95% of the ground material passes through a 20 mesh sieve (about the particle size of fine espresso).

1. If not using Neogen's prepared solution, prepare a 65% ethanol solution by mixing 6.5 parts ethanol with 3.5 parts distilled or deionized water for each sample.
2. Combine 10 grams of ground sample with 20 mL of 65% ethanol. Vigorously shake, using hand or mechanical means, for 1 minute.
3. Allow the sample to settle, then filter with a filter syringe or Whatman No. 4 filter paper.
4. You may also pipette 1 mL of sample into a 1.5 mL micro-centrifuge tube, and centrifuge for 1 minute using a microcentrifuge.
5. The sample is now ready for testing.

TEST PROCEDURE

1. Place the appropriate number of red sample cups into a sample cup rack.
2. Add 100 μ L of the clear portion of the sample extract (or filtered extract) to the red sample cup.
3. Using a new tip, add 500 μ L of sample diluent to the red cup containing the sample extract. Mix well by pipetting up and down 3 times.
4. Place the Reveal for Aflatoxin test strip into the red sample cup and set a timer for 3 minutes. Ensure the test strip comes into contact with liquid and begins to wick.
5. Remove the strip from the sample cup after it has developed for 3 minutes.
6. Read visually by holding strip up to line of sight. Strip must be read within 30 seconds of the 3 minute incubation.

INTERPRETATION OF RESULTS

Results less than 20 ppb: If a line forms in the test zone and another line forms in the control zone within 3 minutes (2 lines total), the sample contains less than 20 ppb aflatoxin (negative sample). A negative sample may be determined as soon as 2 lines are visible on the strip. If there is no line in the control zone, the test is invalid and the sample should be re-tested with another strip.

Results greater than or equal to 20 ppb: If after the full 3 minutes there is no visible line in the test zone, but a visible line in the control zone, the sample contains at least 20 ppb aflatoxin (positive sample). If there is no line in the control zone, the test is invalid and the sample should be retested with another strip.


Notes:

- Observations should be made between 3 and 3.5 minutes. Observations after 3.5 minutes may be inaccurate due to over-development of the device.
- The strips can be read using Neogen's Reveal AccuScan Gold Reader*. This eliminates subjectivity by interpreting and storing sample results.
- Positives can be confirmed by Neogen's quantitative Veratox* products for aflatoxin or HPLC.

*Contact Neogen or visit www.neogen.com for more information on these products.

PERFORMANCE CHARACTERISTICS

Reveal for Aflatoxin is designed to screen for aflatoxin in corn at 20 ppb.

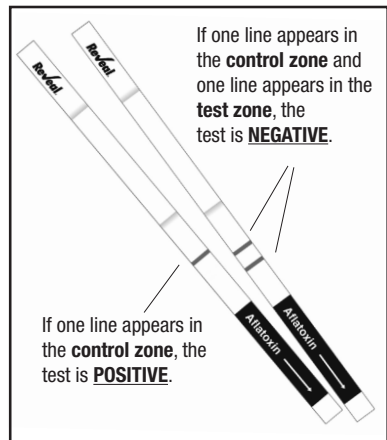Neogen Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all Neogen test kits, is available.

SDS INFORMATION AVAILABLE

Safety data sheets (SDS) are available for this test kit, and all of Neogen's test kits, on Neogen's website at foodsafety.neogen.com, or by calling Neogen at 800/234-5333 or 517/372-9200.

TERMS AND CONDITIONS

For Neogen's full terms and conditions, please visit www.neogen.com/en/terms-and-conditions.

WARRANTY

Neogen Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, Neogen will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. Neogen shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

TEST KITS AVAILABLE FROM NEOGEN

Natural toxins

- Aflatoxin, DON, ochratoxin, zearalenone, T-2/HT-2 toxins, fumonisin, histamine

Foodborne bacteria

- *E. coli* O157:H7, *Salmonella*, *Listeria*, *Listeria monocytogenes*, *Campylobacter*, *Staphylococcus aureus*, *Salmonella enteritidis*

Sanitation

- ATP, yeast and mold, total plate count, generic *E. coli* and total coliforms, protein residues

Food allergens

- Almonds, crustaceans, eggs, gliadin, hazelnut, milk, mustard, peanuts, sesame, soy, walnuts, multi-treenut

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