

Read instructions carefully before starting test

Reveal[®]

for Horse

For the qualitative detection of raw horse meat

REFRIGERATE AT 2–8°C (35–46°F) • DO NOT FREEZE

SPECIES IDENTIFICATION

Preventing adulteration of feed and food produce with non desirable or low quality animal products is important for economic, regulatory, health, and ethnic reasons. The identification of species is performed to assure consumers that the products they purchase are safe, wholesome, and properly labelled.

INTENDED USE

Reveal for Horse is an immunochromatographic lateral flow assay used for the rapid, qualitative detection of horse meat in raw meats and processed raw meat products, such as sausages or burgers. The test can detect as little as 0.5% horse meat in a non-target meat matrix.

Reveal[®] for Horse is intended for use by quality control personnel or other personnel familiar with handling meat samples only in an industrial food manufacturing or preparation context. Because of the problems of adequate sampling and extraction of meats, it is not suitable for the testing of foods to be consumed in the home or in a restaurant by consumers.

LIMIT OF DETECTION

The test is highly sensitive and specifically designed to screen for low levels of raw horse in raw meat or raw meat products. The limit of detection of the test is 0.5% raw horse meat in a non-target raw meat matrix.

Using environmental swabs, levels of 150 µg/100 cm² of horse protein can be detected.

When analyzing rinses, horse protein residues are detectable at a level of 10 µg/mL. The presence of cleaners and sanitizers can affect limit of detection (LOD) in rinses.

Reveal for Horse has been tested on a panel of raw meat cross-reactants. Cross-reactivity is expected in raw meat from the Equidae family, comprising of horses, donkeys, and zebra. No cross-reactivity against meat from non-target species such as beef, pork, poultry, or sheep was observed. Reveal for Horse validation report is available on request, please contact your NEOGEN[®] representative.

ASSAY PRINCIPLES

Reveal® for Horse is a single step lateral flow immunochromatographic assay. The sample extract is wicked through a reagent zone, which contains antibodies specific for horse serum albumin proteins conjugated to coloured particles. If horse serum albumins are present, they will be captured by the conjugated antibodies. The horse antibody particle complex is then wicked onto a membrane which contains a zone of antibody specific for horse proteins. This zone captures the complex allowing the particles to concentrate and form a visible line. If no horse is present, no line will form. The membrane also contains a control zone where an immune complex present in the reagent zone is captured by an antibody, forming a visible line. The control line will always form regardless of the presence of horse, ensuring the strip is working properly.

MATERIALS PROVIDED

1. 10 Reveal for Horse lateral flow test strips
2. 10 clear sample cups
3. 10 vials of sample diluent
4. 10 disposable 100 µL fixed volume pipettes
5. 10 disposable 200 µL fixed volume pipettes

MATERIALS REQUIRED BUT NOT PROVIDED

1. Extraction containers to contain 10 mL deionized or distilled water for extraction
2. Deionized or distilled water
3. Reveal sample cup rack (NEOGEN® item 9475)
4. Timer (NEOGEN item 9426)
5. Mincer or knife (to finely chop sample)
6. Scale capable of weighing 0.5–400 g ± 0.1 g (NEOGEN item 9427)
7. Measuring cylinder capable of measuring 10 mL
8. Reveal for Meat Speciation Swabbing Kit (NEOGEN item 9541, contains 50 swabs)

STORAGE

Store kit components refrigerated at 2–8°C (35–46°F) to ensure full shelf life. Do not freeze. Test strips should remain capped in their original sample tubes until used to ensure optimal performance.

LIMITATIONS

A negative test cannot exclude the possibility that the food contains horse since it may be distributed unevenly in the food product and may be below the detection limit of the test with that specific sample.

Reveal for Horse is a qualitative test and should only be used as a preliminary screen for the presence of raw horse. This test is not suitable for cooked meats. Due to variations in processing, cured and fermented meat products should be validated prior to routine use. The validity of results obtained with the test should preferably be viewed in conjunction with data from an alternative method, such as an enzyme-linked immunosorbent assay (ELISA) or polymerase chain reaction (PCR).

PRECAUTIONS

1. The test strips must remain inside the stay dry tube before use.
2. Store test kit refrigerated at 2–8°C (35–46°F) when not in use. Do not freeze.
3. Remove the test kits and allow to equilibrate at room temperature (18–30°C, 64–86°F) before use (20–30 minutes out of refrigerator).
4. Do not use kit contents beyond expiration date.
5. Gloves and other protective apparel should be worn at all times to handle raw meat and during test procedure.
6. To avoid cross-contamination, use clean pipettors, extraction container, and sample diluent vial for each sample.

EXTRACTION PROCEDURE FOR RAW MEAT SAMPLES

The sample should be collected according to accepted sampling techniques.

1. Obtain a representative meat sample. Mince or finely chop meat sample until homogenous.
2. Add 1 g of prepared meat sample to 10 mL of distilled or deionized water.
Note: Following validation, the extraction of larger sample sizes may be applicable when using an extraction ratio of 1 part meat sample to 10 parts distilled or deionized water.
3. Vigorously shake for 1 minute.
4. Use the sample supernatant as the sample extract for testing.
5. Discard extracts after completion of analysis.


TEST PROCEDURE FOR RAW MEAT SAMPLES

1. Take the appropriate number of sample diluent vials and sample cups from the test kit. Label cups and vials on the outside for identification if necessary.
2. Remove the cap from the sample diluent vial. Using an 200 μ L fixed volume pipette, transfer 200 μ L of the sample supernatant into the sample diluent vial.
3. Re-cap the sample vial and mix by shaking for 10 seconds.
4. Using an 100 μ L fixed volume pipette, transfer 100 μ L of the sample extract from the sample diluent vial into a new clear sample cup.
5. Place a new Reveal[®] for Horse test strip with the sample end down into the sample cup and set a timer for 5 minutes. Ensure the test strip comes into contact with the liquid and begins to wick.


EXTRACTION AND TEST PROCEDURE FOR SWABS

Note: Swabs are not included in the test kit and can be ordered separately (NEOGEN[®] item 9541).

1. Estimate a swabbing area of approximately 10 cm x 10 cm. Alternatively, use the swab to collect samples of contamination from problem areas (e.g., of processing equipment).

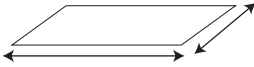


2. Gather the sample with the swab, using one of the following methods:
For dry surfaces: Remove a sterile swab from the packaging and wet the surface with 100 μ L deionized or distilled water using a disposable 100 μ L fixed volume pipette. Swab a 10 x 10 cm area by using a crosshatch technique revolving the swab on the surface. Repeat this swabbing procedure using movements at right angles to those used in the first swabbing.
For wet surfaces: Remove a sterile swab from the packaging and swab a 10 x 10 cm area by using a crosshatch technique revolving the swab on the surface. Repeat this swabbing procedure using movements at right angles to those used in the first swabbing. Do not moisten surface prior to use.

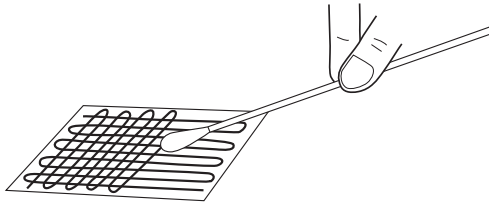

3. Place the swab into the sample diluent vial and carefully break off the moistened end at the prescored mark so that it remains in the vial.

- Secure the cap of the sample diluent vial, to ensure that the stem does not prevent the tube from being properly sealed. Shake for 1 minute.
- Using a disposable 100 μL fixed volume pipette, transfer 100 μL of the sample extract from the sample diluent vial into a new sample cup.
- Place a new Reveal for Horse test strip with the sample end down into the sample cup and set a timer for 5 minutes. Ensure the test strip comes into contact with the liquid and begins to wick.

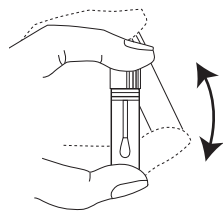

EXTRACTION AND TEST PROCEDURE FOR RINSE WATERS

- Obtain a representative rinse water sample.
- Using a 200 μL disposable fixed volume pipette, add 200 μL of the rinse water sample and transfer into a sample diluent vial.
- Secure the cap of the sample diluent vial and shake for 10 seconds.
- Using a disposable 100 μL fixed volume pipette, transfer 100 μL of the sample extract from the sample diluent vial into a new sample cup. Place a new Reveal[®] for Horse test strip with the sample end down into the sample cup and set a timer for 5 minutes. Ensure the test strip comes into contact with liquid and begins to wick.
- Place a new Reveal for Horse test strip with the sample end down into the sample cup and set a timer for 5 minutes. Ensure the test strip comes into contact with the liquid and begins to wick.

READING RESULTS

Test results should be read within 1 minute of completion of the 5 minute incubation.

Negative Result


No line at the test line position: Level of horse is below detection limit (see limitations section) or not present in the sample. Control line present.

Positive result

Any intensity of a red line at the test line position: Level of horse is above detection limit. Control line present.

Invalid results

If a red line does not appear at the control line position, then the test is invalid.

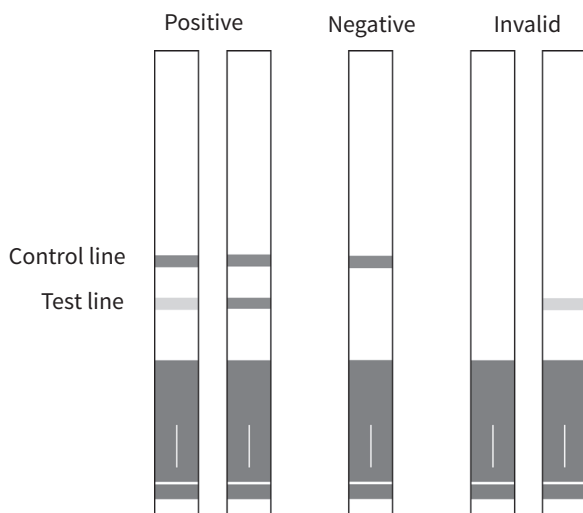

CUSTOMER SERVICE

NEOGEN® Customer Service and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all NEOGEN test kits, is available.

SDS INFORMATION AVAILABLE

Safety data sheets (SDS) are available for this test kit, and all of NEOGEN's test kits, on NEOGEN's website at neogen.com, or by calling NEOGEN at 800.234.5333 or 517.372.9200.

TERMS AND CONDITIONS

For NEOGEN's full terms and conditions, please visit neogen.com/en/terms-and-conditions.

WARRANTY

NEOGEN Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, NEOGEN will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. NEOGEN shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

TESTING KITS AVAILABLE FROM NEOGEN

Natural Toxins

- Aflatoxin, ergot alkaloids, DON, ochratoxin, zearalenone, T-2/HT-2 toxins, fumonisin, histamine

Foodborne Bacteria

- *E. coli* O157:H7, *Salmonella*, *Listeria*, *Listeria monocytogenes*, *Campylobacter*, *Staphylococcus aureus*, *Salmonella enteritidis*

Sanitation

- ATP, yeast and mold, total plate count, generic *E. coli* and total coliforms, protein residues

Food Allergens

- Almonds, coconut, crustaceans, eggs, gliadin, hazelnut, milk, mustard, peanuts, sesame, soy, walnuts, multi-treenut

Genetic Modification

- CP4 (Roundup Ready®)

Ruminant By-products

- Meat and bone meal, feed

Species Identification

- Raw and cooked meat samples



North America

NEOGEN Headquarters

800.234.5333 (U.S./Canada)
foodsafety@neogen.com
NEOGEN.com

Europe, Middle East, and Africa

NEOGEN Europe

+ 44 (0) 1292 525 600
info_uk@neogeneurope.com
NEOGEN.com

Mexico

NEOGEN Latinoamerica

+52 (55) 5254-8235
informacion@neogenlac.com
NEOGEN.com

Brazil

NEOGEN do Brasil

+55 19 3935.3727
info@neogendobrasil.com.br
NEOGEN.com

China

NEOGEN Bio-Scientific Technology

+86 21 6271 7013
info@neogenchina.com.cn
www.neogenchina.com.cn

India

NEOGEN Food and Animal Security

+91 484 2306598, 2301582
info@neogenindia.com
www.neogenindia.com

©NEOGEN Corporation, 2020. NEOGEN and Reveal are registered trademarks of NEOGEN Corporation. All other brand and product names are trademarks or registered trademarks of their respective companies.

Patent: <https://www.neogen.com/patents/>

9531B

Official_9531B_9531_Reveal_Horse_ES