

Reveal[®] 3-D for Gliadin

DOWNLOAD AND READ KIT INSTRUCTIONS COMPLETELY BEFORE PERFORMING TEST.



Materials Provided:

- 20 Reveal[®] 3-D for Gliadin test devices
- 20 bottles of extraction buffer
- 20 sterile swabs
- 01 bottle swab wetting solution

Product Number: 8519

- Rinse: 5 ppm
- Swab: 1 µg/100 cm²
- Testing time: 5 minutes

Call 800.234.5333 to order or visit neogen.com

Rinse Sampling




Add 0.25 mL of the sample to an extraction buffer bottle.



Secure cap and shake for 1 minute. Proceed to swab/rinse testing step 1.

Swab Sampling

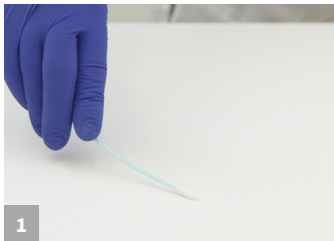

Gather the sample. For dry surfaces, moisten with swab wetting solution. Do not premoisten for wet surfaces.



Insert the swab in the extraction buffer bottle and break off into the bottle.



Secure cap and shake for 1 minute. Proceed to swab/rinse testing step 1.

Swab/Rinse Testing



Remove the lid and fill with liquid from the bottle.



Dip the Reveal[®] 3-D device into the liquid within the lid, ensuring the liquid saturates the cavity.



When the liquid is seen running up the test window, place the device on a flat surface. Allow the test to develop for 5 minutes.



Interpret the results.

